

**PHOSPHATE BUFFERED SALINE (PBS)****PHOSPHATE BUFFERED SALINE (PBS)**

Concentrated PBS (10x)

INbusal

Expiration date > 8 years
Store at room temperature

**SUMMARY****REAGENTS**

INbusal Concentrate, product number 2134, is liquid. Store at room temperature.

Sodium Chloride

Disodium phosphate

Monosodium phosphate

For stabilities refer to expiry date printed on the product label.

PROCEDURE

Dissolve the contents of one vial INbusal (2134) with 9.0 l distilled or deionised water in a INbusal container (2029) and mix carefully. For smaller wished amounts of PBS dilute and mix 1 part INbusal concentrate (2134) with 9 parts distilled or deionised water. The pH will be 7.40.

PRODUCTS

Products	Product no.	Quantity
INbusal Concentrate	2134	1 x 1 liter
INbusal Container	2029	1 x 10 liter

**NOTES**

1. For in vitro diagnostic use only.
2. For professional use only.
3. Contact INstru chemie for the latest edition product insert.
4. Printed in the Netherlands, July 2015, version 4.0