



INSTRUCLEAR - WATERCLEAR

ANTISEPTIC CONCENTRATE TO MAINTAIN THE QUALITY OF WATER IN THERMOSTATS, WATER BATHS, CUVETTES SYSTEMS, ETC.

INstruclear – Waterclear:
Visual blue colour indicator
Ready for use
Stability > 6 years after production



SUMMARY

REAGENTS

INstruclear, product number 2316, is an antiseptic concentrate. Store at room temperature.

PRINCIPLE

INstruclear keeps water germ free and is complex binding active, INstruclear prevents grow of algae, moulds and bacteria, bad smell, calcium and heavy metal formations.

Protection against stoppage and pollution of equipments, tubes, cuvettes, etc. and prevention of the users against mould and/or bacterial infections. The optimal conduction of heat is guaranteed during several years.

INstruclear saves work and costs because the water only has to be renewed if the water stays turbid. INstruclear protects your instruments and instrument tubing against dirt, blocking and strong wear and tear. INstruclear helps the constant conduction of heat by keeping the instruments parts clear.

PROCEDURE

Add 5 – 10 ml INstruclear per liter distilled or deionized water into a previous cleaned water bath, thermostat and or relevant equipments.

When the blue colour disappears add again 5 – 10 ml INstruclear per liter into the water bath. The water should be clear and blue. If the water stays turbid after adding a few doses of INstruclear clean the water bath and renew the distilled or deionized water and INstruclear.

PRODUCTS

Products	Product no.	Quantity
INstruclear	2316	6 x 100 ml

NOTES

1. For in vitro diagnostic use only.
2. For professional use only.
3. Always contact INstru**chemie** for the complete product insert and latest edition.
4. Printed in the Netherlands, July 2015 – version 3.0

